DISCOVERY

PRO

ENGLISH	1 – 14
DEUTSCH	15 – 30
FRANÇAIS	31 – 46
ESPAÑOL	47 – 62
PORTUGUÊS	63 – 76
POLSKI	77 – 92
РУССКИЙ	93 – 108

CE IVD

HTL
HTL LAB SOLUTIONS

PZ HTL S.A. is in possession of a registered and certified quality management system which includes the development, production, and sales of high quality Liquid Handling products.

CONTENTS

1. INTRODUCTION
1.1 PRODUCT DESCRIPTION
1.2 WORK SAFETY INSTRUCTIONS
1.3 TECHNICAL PARAMETERS
2. OPERATION
2.1 SETTING THE VOLUME
2.2 PREPARATION FOR OPERATION
3. PIPETTING RECOMMENDATIONS
4. ASPIRATING AND DISPENSING LIQUIDS
4.1 ASPIRATING LIQUIDS
4.2 DISPENSING LIQUIDS
4.3 PRE-RINSING
4.4 ASPIRATING HIGH DENSITY LIQUIDS
5. RECALIBRATION
6. MAINTENANCE
7. TROUBLESHOOTING
8. PIPETTE KIT

1. INTRODUCTION

DISCOVERY PRO pipettes are designed for general laboratory use. Pipettes can be used for measurement and transfer of aqueous solution, acids, bases and enzyme assay applications.

These variable volume pipettes cover the range from 1 μ l to 300 μ l in 8 models.

DISCOVERY PRO pipettes operate utilizing the air cushion principle, i.e. the aspirated liquid has no contact with the shaft or plunger of the pipette. Liquid is aspirated into disposable tips put on the pipette.

Pipettes are equipped with a four-digit counter displaying the set volume, and the aspirated volume is set by means of the adjustment knob (fig. 1B). The pipette design allows the user to lock the volume setting.

1.1. PRODUCT DESCRIPTION

A. Pipetting pushbutton

B. Volume adjustment knob

Apart from adjusting the volume, the volume adjustment knob is also designed to lock the volume setting.

C. Ejector button

D. Cap

It is used to cover the calibration adjustment mechanism and it can be used for pipette identification.

E. Calibration switch

It is used to switch the pipette from operation mode to calibration mode.

F. Volume counter

The four-digit volume counter displays the volume setting.

G. Counter cap

The cap has a printed range of aspirated volume, appropriate for the given pipette model.

H. Tip ejector

I. Shafts

Made of high quality material provides high chemical and mechanical strength.

J. Module

The module consists of pistons assembly and shafts assembly which move independently.

Pipette volume is identified by the color of the calibration switch (fig. 1E) and the cap (fig. 1D), depending on the tip type.

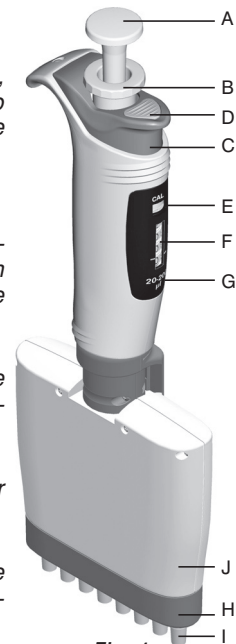
Fig. 1

The applied colors:

- 10 µl pipettes – red,
- 50 µl, 200 µl pipettes – yellow
- 300 µl pipettes - green

1.2. WORK SAFETY INSTRUCTIONS

Used symbols:

 **WARNING!** Risk of injury

CAUTION: Potential damage to the device or pipetting errors

Long-lasting and faultless working of the pipette depends to a large extent on its operation. Read the pipette operating instructions carefully and comply with the principles included.

CAUTION:

- Pipette is designed to work with the tips only. Do not aspirate liquids without tips attached. Aspirated liquid should not get into the pipette because this may damage it.
- Use only the pipette with the tips attached.
- One time use of a tip guarantees safety and eliminates possible contamination of the aspirated liquid.
- Keep the pipette clean; do not cleanse it with chemically aggressive substances (e.g. acetone).
- Do not hold the pipette horizontally if there is liquid in the tip.
- Only use pipettes in accordance with the manufacturer instruction to ensure the proper parameters of pipettes.
- In case of improper operation of the pipette, clean the pipette carefully following the instructions given or send it to the service representative.

 **WARNING!**

- Follow the laboratory work safety regulations.
- Exercise extreme caution while pipetting chemically aggressive substances. Use the protective equipment, glasses and gloves.
- Never direct the pipette with the tips attached towards yourself or anyone else when there is liquid in it.
- Use only the spare parts and accessories recommended by the manufacturer.

1.3. TECHNICAL PARAMETERS

DISCOVERY PRO pipettes are high quality laboratory instruments, which provide the highest measurement accuracy and precision.

Accuracy errors and repeatability of liquid volume measurements depend on quality of the applied tips. The errors provided, have been determined with the use of **HTL** tips by means of gravimetric method based on minimum 10 samples of distilled water with temperature of 20°±1°C under the EN ISO 8655 standard.

Model	Cat.no.	Volume [µl]	Accuracy [%]	Precision [%]	Increment [µl]	Fits to tips [µl]
DP8-10 DP12-10	4951 4955	Min 0.5	± 10.0	± 8.0	0.01	10
		5	± 4.0	± 2.0		
		Max 10	± 2.0	± 1.2		
DP8-50 DP12-50	4952 4956	Min 5	± 4.0	± 2.5	0.05	200
		25	± 3.0	± 1.2		
		Max 50	± 1.6	± 0.6		
DP8-200 DP12-200	4953 4957	Min 20	± 3.0	± 1.5	0.2	
		100	± 1.5	± 0.8		
		Max 200	± 1.0	± 0.6		
DP8-300 DP12-300	4954 4958	Min 50	± 1.6	± 1.5	0.5	300
		150	± 1.2	± 1.0		
		Max 300	± 1.0	± 0.6		

Table 1


Only application of **HTL** tips guarantees accuracy of measurements at the level shown in the table above.

The pipette design enables the user to perform calibration process according to the rules presented in section 5.

2. OPERATION

2.1. SETTING THE VOLUME

Setting the volume is performed by the volume adjustment knob (fig. 1B).

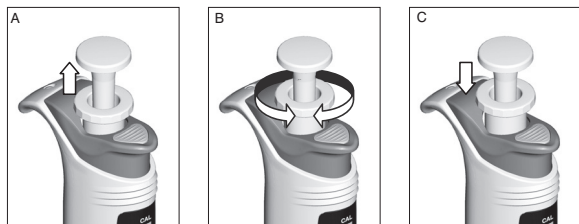
Fig. 2

The range of aspirated volume is displayed on the counter cap (fig. 1G). The volume is displayed by the counter and consists of four digits, which should be read from top to bottom. The digits on the lower drum constitute minimum graduation for the particular model. Value of the minimum graduation (increment) is presented in Table 1.

Volume can be changed after choosing the upper position of the knob (fig. 2A and 2B).

After setting the desired volume, the knob (fig. 1B) should be locked by pushing into the lower position (fig. 2C).

Examples of counter indications: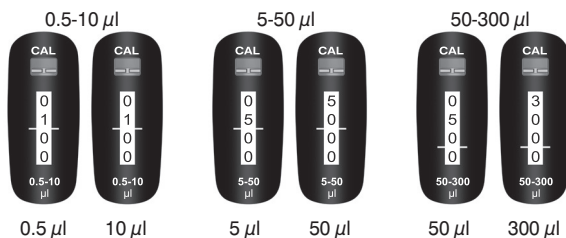


Counter indications above the horizontal indicators shown volume in [µl].

To achieve maximum accuracy, the required volume should be set from higher volume by reducing indications of the counter.

Example of correct volume setting:

- If the required volume is lower than the volume set on the counter it is necessary to reduce indications of the counter to the required value. Before achieving the required value it is necessary to reduce the speed

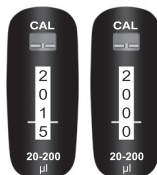Fig. 3

with which you rotate the adjustment knob (fig. 1B), so as not to exceed the value (fig. 3).

- If the required volume is higher than the volume set on the counter it is necessary to increase indications of the counter to the value which exceeds the required volume by rotating the bottom drum by about 1/3 of its circumference. Then, by slowly rotating the adjustment knob (fig. 1B), reduce the setting to the required value (fig. 3).

After setting the required volume, the knob (fig. 1B) should be moved downward to lock the set volume (fig. 2C).

CAUTION: The calibration switch (fig. 1E) which is above the counter should remain in the lower position during operation. Otherwise, each change of capacity will result in decalibration of the pipette.

2.2. PREPARATION FOR OPERATION

Place the tips on the shafts of the pipette. Holding the pipette vertically, press it against the tips in the rack box, until the shafts retreat about 1.5 mm into the multichannel module. The suspension system of the shafts ensures even and tight sealing of the tips.

The twisting motion does not have to be performed.

See Table 1 for the appropriate tips designed for the given pipette model.

CAUTION:

- It is recommended to use the tips recommended by the manufacturer, as only those tips will ensure accurate and repeatable liquid measurements.
- Never aspirate liquid into a pipette without tips attached.

3. PIPETTING RECOMMENDATIONS

Observing the following recommendations will ensure accuracy and repeatability of liquid sampling.

- While operating **DISCOVERY PRO** pipette the volume setting should be locked - the adjustment knob in its lower position (fig. 2C).
- Operate the pipette pushbutton slowly and smoothly when liquid aspirating and dispensing.
- Pipette is equipped with a switch of counter's mode of operation – calibration switch (fig. 1E). This switch should be shifted only for the time of calibrating the pipette (upper position marked with the "CAL" symbol).

- The depth of tips immersion in the aspirated liquid should be around 2-4 mm.
- While operating, the pipette should be held in a vertical position.
- New tips should be pre-rinsed.
- The tips should be replaced with new ones if:
 - a different type of liquid is handled;
 - a different liquid volume is dispensed;
 - visible droplets of liquid remain in them.
- Pipetted liquid should not enter the pipette shafts. To ensure this:
 - press and release the pipetting pushbutton slowly and smoothly;
 - do not put the pipette away if there is any liquid in the tips;
 - do not invert the pipette;
 - do not set and aspirate volumes exceeding nominal values.
- Before aspirating liquid of temperature that differs from ambient temperature, it is recommended to rinse the tips several times with the aspirated liquid.

4. ASPIRATING AND DISPENSING LIQUIDS

4.1 ASPIRATING LIQUIDS

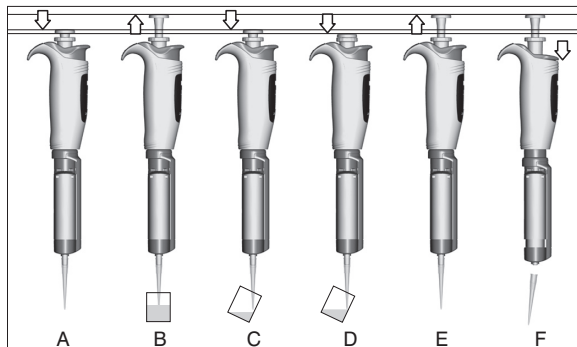
Fig. 4

- The pipetting pushbutton (fig. 1A) should be pressed to the first stop (fig. 4A),

- While holding the pipette in a vertical position, tips should be immersed into the aspirated liquid to the depth of 2 to 4 mm, and flushed once by drawing a dose of liquid and dispensing it out with slow and steady movement. The tips may aspirate air if immersed to a depth lower than recommended or if the pipetting pushbutton is released quickly,
- To aspirate liquid release the pipetting pushbutton smoothly (fig. 4B),
- Wait for about 1 second before withdrawing the tips from the aspirated sample, fig. 4.

⚠ WARNING! Do not touch the used tips.

4.2 DISPENSING LIQUIDS

- While holding the pipette slightly inclined from 10 to 40° off the vertical, place the tips orifice against the inside wall of the receiving vessel.
- Then, press the pipetting pushbutton smoothly to the first stop, dispensing the liquid (fig. 4C).
- After about one second press the pipetting pushbutton to the final stop, blowing out the remaining liquid from the tips (fig. 4D).
- Holding the pushbutton depressed remove the pipette, drawing it against the vessel inner wall.
- Release the pipetting pushbutton to its starting position (fig. 4E) and eject the tips by pressing the tip ejector button (fig. 4F).

CAUTION: Replace the tips with a new one whenever a different liquid is to be pipetted.

4.3 PRE-RINSING

When pipetting liquids of viscosity higher or surface tension lower than those of water (e.g. serum or organic solvents) a film of liquid is formed on the inside wall of the pipetting tips. The film can create an error. Since the film remains relatively constant in successive pipetting operations with the same tips, this error can be avoided by forming the film before the first pipetting. This is done by performing a full pipetting cycle of a liquid into the same vessel. After such a procedure, the film of liquid is already formed in the tips and ensures better accuracy and repeatability of successive pipetting operations.

4.4 ASPIRATING HIGH DENSITY LIQUIDS

The accuracy errors provided in Table 1 have been determined using distilled water. When pipetting liquids whose properties (density, viscosity, surface tension) differ from the properties of water, it may be necessary to compensate for volume settings.

CAUTION: Aspiration differences can be avoided if the pipetting operation is conducted slowly, so that the liquid can slowly adjust to the changing pressure. Wait for about 2 seconds after each aspirating and dispensing, without changing the position of the pipette.

If the above procedure does not provide accurate results, pipette recalibration is required according to section 5.

After calibration, it is recommended to record recalibration and the adjustment value, which will facilitate return calibration to the standard liquid.

5. RECALIBRATION

DISCOVERY PRO pipettes are calibrated by gravimetric method, using **HTL** tips and distilled water, at the temperature $20^{\circ} \pm 1^{\circ} \text{C}$, according to EN ISO 8655 standard.

It is recommended to conduct periodic inspection of the pipette operation, with the interval between inspections being dependent on the type of load (aspirated liquids) and other conditions (load intensity and autoclave sterilization, replacement of subassemblies) in which the pipette is used.

The pipette recalibration procedure should be carried out if during the pipette operation you find that the accuracy error (the difference between the real aspirated volume and the preset volume) exceeds the permissible value given in Table 1.

Aspiration efficiency of the pipette is determined e.g. by: applied tips, properties of pipetted liquid (density, viscosity) and operation conditions (ambient temperature, pressure).

Checking the pipette parameters

To determine accuracy error of the pipette, the following requirements must be fulfilled:

- The ambient temperature of the pipette, tips and the test liquid (typically distilled water) should be within the range of $20\text{-}25^{\circ}\text{C}$ and stabilized during weighing within the range $\pm 0.5^{\circ}\text{C}$,

- The balance sensitivity should be appropriate to the tested volume (Table 2),
- Table 3 shows the values of the $[\text{mg}/\mu\text{l}]$ conversion rates for distilled water or another liquid of comparable density.

Volume $[\mu\text{l}]$	0.1 - 10	10 - 100	> 100
Balance sensitivity $[\text{mg}]$	≤ 0.001	≤ 0.01	≤ 0.1

Table 2

Temperature $^{\circ}\text{C}$	Pressure $[\text{kPa}]$		
	95	101.3	105
20	1.0028	1.0029	1.0029
21	1.0030	1.0031	1.0031
22	1.0032	1.0033	1.0033
23	1.0034	1.0035	1.0036
24	1.0037	1.0038	1.0038
25	1.0039	1.0040	1.0040

Table 3

CAUTION: The pipetting procedure should be conducted in accordance with the rules described in sections 3 and 4.

Recalibration procedure:

- Set the dose volume depending on the pipetted volume according to Table 4.

Model	Range of the pipettes volumes $[\mu\text{l}]$	Preset volume $[\mu\text{l}]$	Permissible volumes $[\mu\text{l}]$
DP8-10, DP12-10	0.5-10	0.5	0.45-0.55
DP8-50, DP12-50	5-50	5	4.8-5.2
DP8-200, DP12-200	20-200	20	19.4-20.6
DP8-300, DP12-300	50-300	50	49.2-50.8

Table 4

- Perform 5 aspirations, weigh each one and calculate the average value of the aspirations.
- Calculate average aspirated volume in $[\mu\text{l}]$ by multiplying the average aspiration amount $[\text{mg}]$ by the distilled water density coefficient $[\mu\text{l}/\text{mg}]$, which depends on temperature and pressure (Table 3).
- If the average aspirated volume exceeds the permissible values, pipette recalibration should be conducted.

Pipette recalibration consists in setting the real value obtained during weighing on the counter. To do so, the following should be conducted:

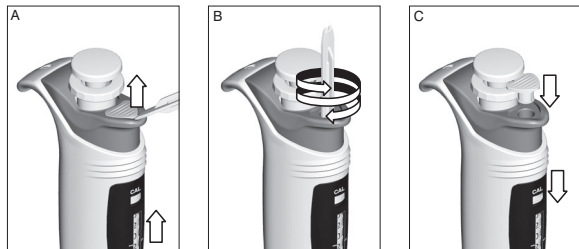
Fig. 5

- Remove the cap (fig. 1D) using the calibration key (fig. 5A),
- Set the calibration switch (fig. 1E) in its upper position (fig. 5A),
- Insert the calibration key into the ejector orifice (fig. 1C) and place it in the calibration screw of the counter (fig. 5B),
- Turn the key so as the volume indicated by the counter is equal to the average calculated volume.
- Remove the calibration key and set the calibration switch in its lower position (fig. 5C), thus going to the pipetting mode,
- Place the cap of the ejector pushbutton (fig. 5C).

After conducted the calibration procedure it is recommended to record recalibration.

CAUTION: Before calibration, the volume adjustment knob should be blocked. If the knob is shifted after switching the pipette on the calibration mode ("CAL"), calibration of the pipette will be changed without a visible change in settings of the counter.

6. MAINTENANCE

The pipette requires periodic maintenance, depending on the applications performed and intensity of use. The subassemblies subject to action of aggressive solution vapors, such as shafts elements, should be regularly inspected and cleaned.

External surfaces of the handgrip, knob and pushbuttons may be cleaned with a tissue dipped in isopropyl alcohol.

CAUTION: Before using cleaning agents other than those recommended by the manufacturer, check the compatibility charts and consider chemical resistance of the following plastics: PP, PC, POM, PA, PPS, PVDF used to produce the pipette parts.

The pipette can be sterilized in the autoclave at the temperature of 121°C for 20 minutes. After sterilization, the pipette should be dried and cooled down to room temperature. It is recommended to sterilize the pipettes in autoclave with initial vacuum and drying cycle.

CAUTION:

- Sterilization in other conditions may cause the damage of the pipette,
- Set the volume adjustment knob in its upper (unlocked) position before sterilization.

Repeatability of obtained results does not change if the pipette is operated properly and the autoclaving procedure is conducted appropriately. As there can be a slight change in the dosing accuracy it is recommended:

- To check calibration of the pipette after 1st, 3rd and 5th autoclaving procedures, and every 10 autoclaving cycles during further operation of the pipette.

The pipettes are UV resistant which was confirmed by our tests. The recommended distance from the radiation source to exposed element should be not less than 50 cm. Prolonged and very intense UV exposure can cause de-coloration of pipette parts, without affecting its performance.

7. TROUBLESHOOTING

If you notice an improper pipette operation, identify the cause and eliminate the fault. To do this, follow the instruction in the sequence provided. Replacement of parts should be required only occasionally, and should not occur under normal pipette use.

Problem	Cause	Instructions
Droplets of liquid remain in the pipette tips	The liquid is dispensed too fast	Decrease the dispensing speed
	The tips wettability has increased due to extensive use	Replace the tips with new ones
Air bubbles appear in the liquid aspirated into the tips	The pipette tips immersion is too shallow	Immerse the tips deeper according to the instruction
	The pipette tips are incorrectly pressed onto the pipette shafts	Press the tips firmly
	The tip is damaged or worn out due to extensive use	Replace the tips with new ones
The pipette incorrectly aspirates the liquid or the liquid drops out from the tips	The pipette tips are incorrectly pressed onto the pipette shafts	Press the tips firmly
	The shafts surface are damaged or contaminated at the sealing site	Clean the shafts or send the pipette for service in order to replace the shafts
	The plungers or the seals are damaged due to prolonged aspiration of aggressive liquids	Send the pipette for service
	The inside of the pipette is contaminated	
	No lubrication on the sealing elements	

Problem	Cause	Instructions
Uneven work of the pipetting set, the pipetting pushbutton gets blocked	The inside of the pipette is contaminated due to aspiration of aggressive substances	Send the pipette for service
	The inside of the pipette is contaminated due to the liquid entering the pipette	
	No lubrication on the sealing elements, e.g. after repeated autoclaving procedures	
Incorrect aspiration	Liquid with properties other than water (density, viscosity)	Calibrate the pipette using the liquid which is to be pipetted
	The volume adjustment knob used in the "CAL" mode	Calibrate the pipette according to section 5

If the problem continues after carrying out the above steps, contact your pipette supplier to conduct repairs.

CAUTION: Before returning the pipette, please ensure that the pipette is completely free of any aggressive chemical, radioactive, microbiological contamination which could pose a threat during transport and repair.

Clean the pipette as far as it is possible.

8. PIPETTE KIT

The pipettes are delivered in the kits including:

- Pipette
- Short instruction manual
- CD
- Calibration key
- Pipette stand

All rights reserved. Product described in this manual is subject to availability and technical modification. Errors excepted.

PZ HTL S.A. reserves the right to improve, enhance or otherwise modify its products without prior notification.

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© 2015 PZ HTL S.A.